

Procedures Mixing Antibiotics for Dermato Tests.

1. Chloromycetin (^{1,000,000 units} 1 gram bottle) - Add
4.5 ml Sterile H₂O will Rehydrate
to 5.0 ml. for Conc. of 200,000 units/ml

2. Rehydrate 1 vial Mycostatin (^{500,000 units} 500,000 units)
(Nystatin) with 10.00 ml. of
Sterile H₂O to give Conc. of 50,000 units/ml.

3. Add $2\frac{1}{2}$ ml of Chloromycetin
mixture to 10 ml mixture
of Mycostatin for Concentration
of $\frac{40,000 \text{ units of N(M) + Chloromycetin}}{12.5 \text{ ml}}$ for $12\frac{1}{2}$ mls of
 $\frac{500,000}{12.5} = \frac{40,000 \text{ units of N+C}}{\text{ml.}}$

4. Add 0.05 ml of 40,000 N+C to
each 10 ml. Tube of Thioglycollate
for Final Conc. of 200 units M+C
per ml of media

LM-WE

Gail Fraser.
4756

2002
19
83

Credit Report.com